J-Box BTLE ISM Sensor Signal Converter

Instruction Manual

–	Regulatory approval	2	ja	取扱説明書	40
cs	Návod k obsluze	4	ko	사용 매뉴얼	44
da	Brugsanvisning	8	nl	Gebruikershandleiding	48
de	Betriebsanleitung	12	pl	Instrukcja obsługi	52
en	Instruction manual	16	pt	Manual de instruções	56
es	Manual de instrucciones	20	ru	Инструкция по эксплуатации	60
fi	Käyttöohje	24	sv	Bruksanvisning	64
fr	Instructions d'utilisation	28	th	คู่มือคำแนะนำ	68
hu	Használati utasítás	32	zh	说明书	72
it	Istruzioni per l'uso	36			

Regulatory approval

FCC Statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy, and if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

IC Statement

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device.

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Regulatory approval

For regulatory approvals please refer to pages 2 and 3.

1 Introduction

The J-Box BTLE (Bluetooth Low Energy) outputs standardized, simulated nA (for oxygen sensors) and mV (for pH sensors) signals. It is a device for the connection of digital ISM sensors (1-wire pH and RS485 optical DO) to an analog transmitter/biocontroller. The J-Box BTLE offers a Bluetooth 4.0 interface for calibration, diagnostics and data logging purposes via iSense or iSense Mobile.

2 Technical data J-Box BTLE

Measurement parameters	1-wire: pH RS485: optical DO
ISM	Full ISM functionality
Transmission frequency	2.4 GHz, Bluetooth 4.0
Certification	CE
Communication	Bidirectional
Range	1 m
Transmission interval	1 s
Power supply	24 V (min. 375 mA)
Operating temperature	-15 to +50 °C
Relative humidity	5 ... 95 % rH (non-condensing)
Environmental temperature	-20 to +140 °C
4 – 20 mA current range	3.5 to 22 mA (NAMUR)
Protection class	IP67

2.1 Technical specification mV output

pH signal output	-500 ... +500 mV ($R_0 = 17 \text{ k}\Omega$)
Signal	mV non T-compensated
Zero point adjusted to	0 mV @ 25 °C
Slope adjusted to	-59.16 mV/pH @ 25 °C

2.2 Technical specifications nA output

Output range	0 ... 5000 nA ¹
Signal	nA non T-compensated
Zero point/Offset	0 nA @ 25 °C
Air current/slope adjusted to	-70 nA @ 25 °C

¹ The J-Box BTLE can only be used with sensors that output a simulated electrochemical signal (nA, mV). Do not use sensors that output current (mA)!

3 Operating elements

3.1 Connections

I.	+24 VDC plug	Power supply 24 V/0.75 A angled plug (P/N 30 323 961)
II.	T82 plug	O ₂ (nA) to connect to biocontroller
III.	K9/K8S	pH (mV) to connect to biocontroller

- | | | |
|-----|----------------------|--|
| I. | pH sensor connector | Male BNC connector for ISM 1-wire pH sensor |
| II. | oDO sensor connector | Male VarioPin for ISM InPro 6860i oDO sensor |


3.2 LED messages

- | | |
|-------------------------|--|
| Power (green LED on) | J-Box BTLE powered up successfully |
| Connected (blue LED on) | iSense Mobile or iSense have successfully connected via Bluetooth to J-Box BTLE. |

4 Pairing to iSense Mobile

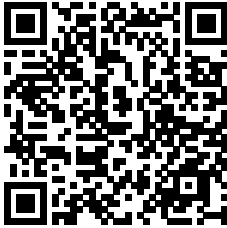The interface works in conjunction with the iSense Mobile app (download from Google Play or Apple App Store) or with iSense PC software.

www.mt.com/isense-mobile

After activating Bluetooth and opening iSense Mobile or iSense, please select the sensor you wish to communicate with.

ISM is a registered trademark of the Mettler-Toledo Group in Switzerland, Brazil, USA, China, European Union, South Korea, Russia and Singapore.

iSense is a trademark of the Mettler-Toledo Group.

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09/2016 © Mettler-Toledo GmbH
France **Mettler-Toledo Analyse Industrielle S.A.S.**
Germany **Mettler-Toledo GmbH**
Switzerland **Mettler-Toledo (Schweiz) GmbH**
Mettler-Toledo GmbH, Process Analytics